

T: +1 (781) 932 3959
F: +1 (781) 932 8705
www.covarisinc.com

Covaris[®]

S220 Series Quick Start Guide

1. Turn on chiller and connect hoses to S220.
NOTE: Ensure proper temperature is selected
2. Fill Covaris tank with de-ionized water to the proper **water level**.
3. Turn on Covaris S220 Series.
4. Turn on laptop.
5. Double Click on the “SonoLab” icon.
NOTE: The software will start and the DEGAS pump will start automatically.
6. Click “NEW” to create a method or the drop-down arrow to load an existing method.
7. Wait 30 minutes for “DEGAS” process to equilibrate and proper temperature to be reached before starting process.
8. Select “RUN” when the system status is ready.
NOTE: Run button will turn green.

COVARIS, INC. 14 GILL STREET UNIT H WOBURN, MASSACHUSETTS, 01801-1721 USA