

PRODUCT INSERT



TITLE: Centrifuge 8 microTUBE Strip V2 Adapter (2)

Check our website for current protocols at <http://covarisinc.com/resources/>



Operational Requirements:

	Name	Part Number
Compatible centrifuge model type	ThermoFisher Scientific mySPIN 12 mini centrifuge	75004081
Centrifuge Rotor Type	Strip Tube Rotor (included with centrifuge)	HS100504
Compatible Covaris microTUBE Strip V2	8 microTUBE-15 AFA Beads Strip V2 (12)	520159
	8 microTUBE-50 AFA Fiber Strip V2 (12)	520174
	8 microTUBE-130 AFA Fiber Strip V2 (12)	520217
Centrifuge Speed	300 RCF maximum	--

Ordering Information:

Name	Part Number
Centrifuge 8 microTUBE Strip V2 Adapter (2)	500541

In this package:

- Centrifuge 8 microTUBE Strip V2 Adapter [Quantity 2]

Revision History:

DCO	Revision	Date	Description of change
021674	A	11/16	Initial Release

Notes:


- Never centrifuge a single 8 microTUBE Strip V2. Always centrifuge two to maintain balance.
- Ensure the 8 microTUBE Strips V2 are fully seated in the adapter prior to spinning.

Installation Instructions:

1. Review the *Rotor Installation and Removal* section of the centrifuge manual. Remove the Strip Tube Rotor (HS100504) and place flat on a bench top.
2. Firmly press the adapter into the middle four holes in the row closest to the finger nut in the center of the rotor. Repeat on the opposite side of the rotor. The adapters should fit tightly and be pressed flush to the bottom.



3. Following the *Rotor Installation and Removal* section of the centrifuge manual, reinstall the rotor into the centrifuge.
4. Load 8 microTUBE Strips V2 into adapters and ensure they are fully seated prior to spinning.

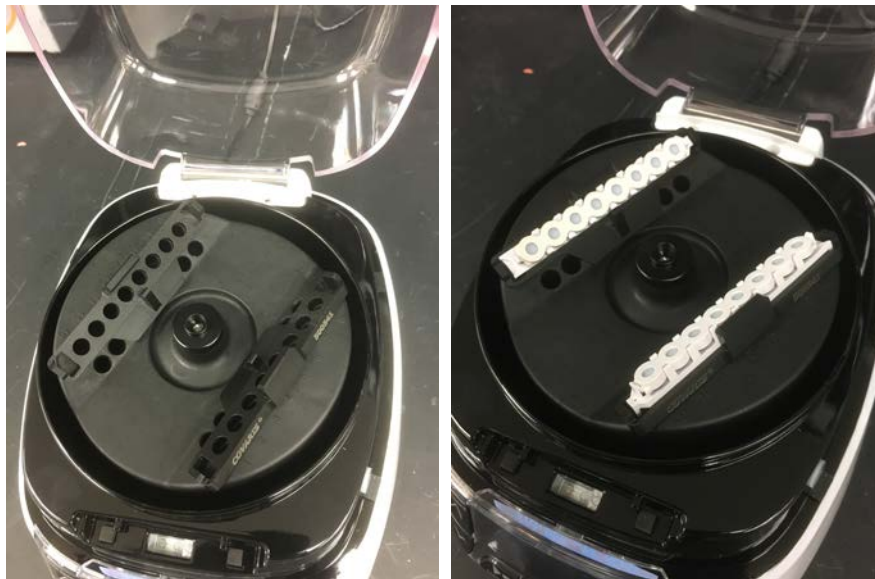

5. Follow instructions provided in manual or protocol for spinning speed and time.

Technical Assistance:

- By telephone (+1 781-932-3959) during the hours of 9:00am to 5:00pm, Monday through Friday, United States Eastern Standard Time (EST) or Greenwich Mean Time (GMT) minus 05:00 hours
- By e-mail at techsupport@covarisinc.com