

PRODUCT INSERT: TT2 Conversion Installation Instructions

Product Warranty Guarantee

Covaris guarantees the performance of all products when used in accordance with our written instruction, under normal operating conditions, and during the expiration period. The user must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any other reason than misuse, Covaris will replace it free of charge. We reserve the right to change, alter, or modify any product to enhance its performance or design. If a product does not meet your expectations, please contact Covaris, Inc.

Technical Assistance

On-going assistance with the operation or application of any of our products is provided via:

- Telephone during the hours of 9AM to 5PM, Monday through Friday, (GMT-05:00) Eastern Time (US & Canada) +1 781 932 3959
- E-mail queries to techsupport@covarisinc.com

INSTRUCTIONS FOR USE

Purpose:

This conversion will allow the Covaris CryoPrep (CP02) to accept larger TissueTubes (TT2).

List of Parts:


1. Cover
2. Tube holder, CryoPrep TT2

List of Tools:

1. 2 mm hex wrench
2. Flat blade screwdriver

Procedure:

1. Replace cover
 - a. Remove tube holder. Center the latch. Close cover.
 - b. Loosen the set screw (1) in the left-rear part of the cover assembly using a 2mm hex wrench. In this picture below, the blue wrench is in position to loosen the set screw.
 - c. Unscrew the cover pivot shaft (2) using a flat-blade screwdriver.
 - d. Install new cover. Reassemble.
2. Install either the TT1 or TT2 tube holder before operating instrument.

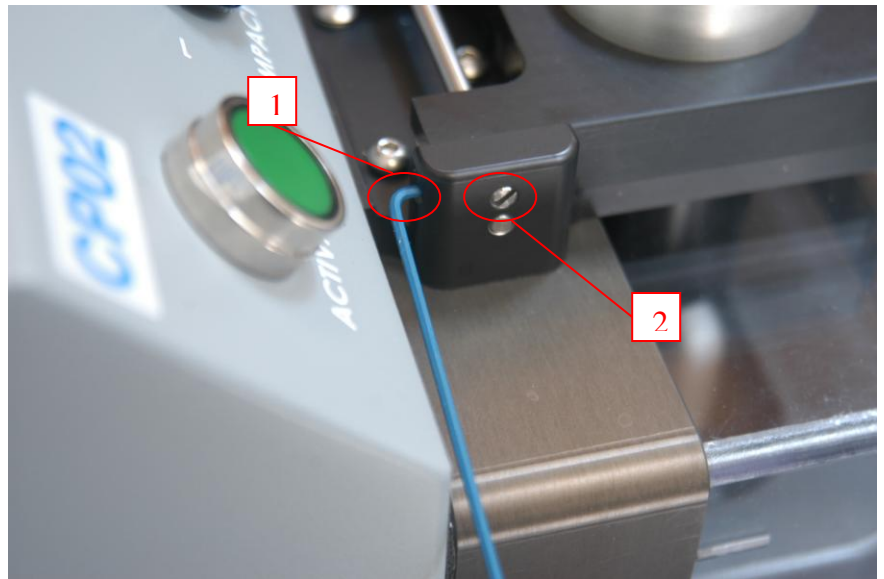

Precaution:



Do not process TT1 tubes in the TT2 tube holder. Replace the TT1 tube holder when processing TT1 parts.