

Procedure for a new install USB drivers and Covaris SonoLab Single™ software on a Windows 7 computer


Note: This install must be performed under an Administrator account.

Under Action Center, click on **change User Account Control Settings**.

Scroll bar to **Never notify**. Click **ok**.

Do you want to allow the following program to make changes to this computer?

Click **Yes**.

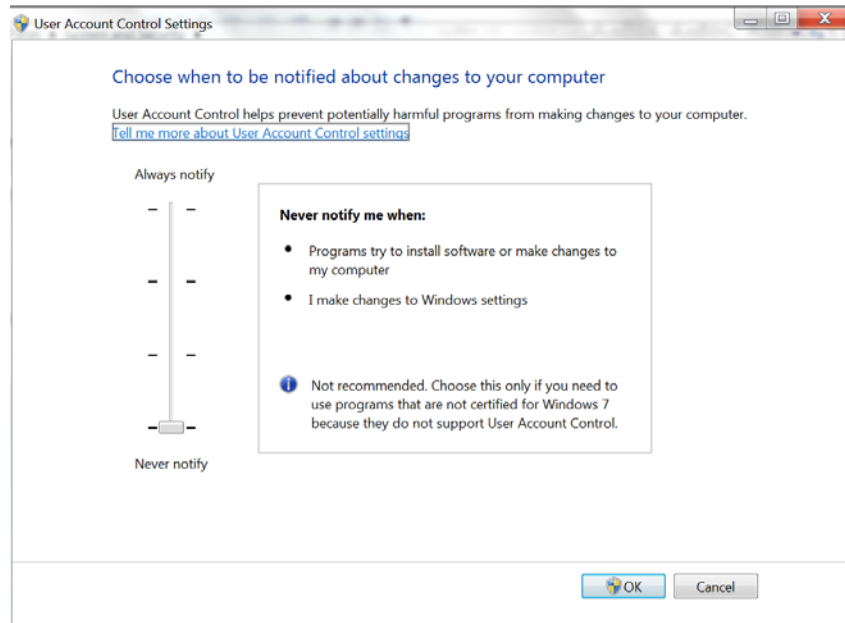

Under Power Options, click on **Change when the computer sleeps**.

In the **Settings for Portable Laptop Power** scheme box select the following for the **On battery** and **Plugged In** section:

Dim the display	Never
Turn Off Display	Never
Put computer to sleep	Never

Save Changes.

Software Installation for the S-Series

Download the appropriate **SonoLab** version and **Drivers** from Covaris and **extract** the Installer and Driver files to the desktop.

Note: At this point turn WIFI OFF.

Right-click “**setup.exe**” in the **SonoLab 2.5.5** folder and **Run as Administrator** to launch the software setup process.

If Windows asks if this is a trusted application, or if you wish to make changes, answer **Yes**.

Follow the instructions for installation. When installation is complete, click on **Finish**.

Hardware Installation for the S-Series

Obtain the instrument and connect it to the computer using a USB Cable and a power cord.


Plug the unit in and power it ON.

Click on the **Start menu\ Control Panel\Hardware and Sound**.


Under **Devices and Printer**, click on **Device Manager**. The Device manager window will appear.

Under “**Other devices**” in the device manager window “**Covaris Sx00 Series**” will appear:


Select **Covaris Sx00**, right click on it and select “**Update Driver Software**”

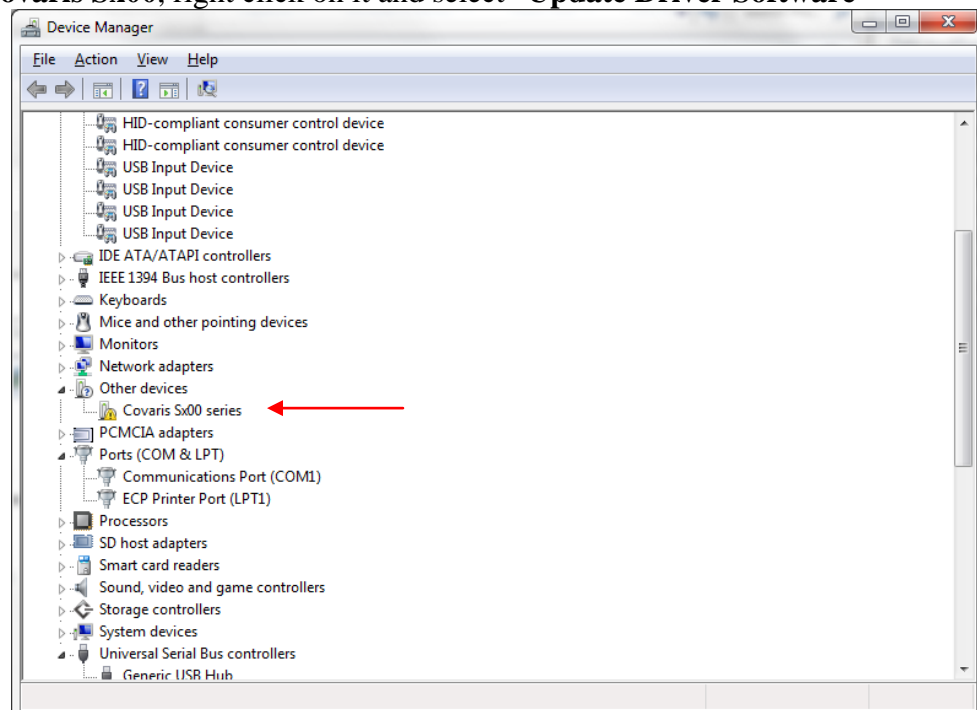


Select **“Browse my computer for driver software”**.

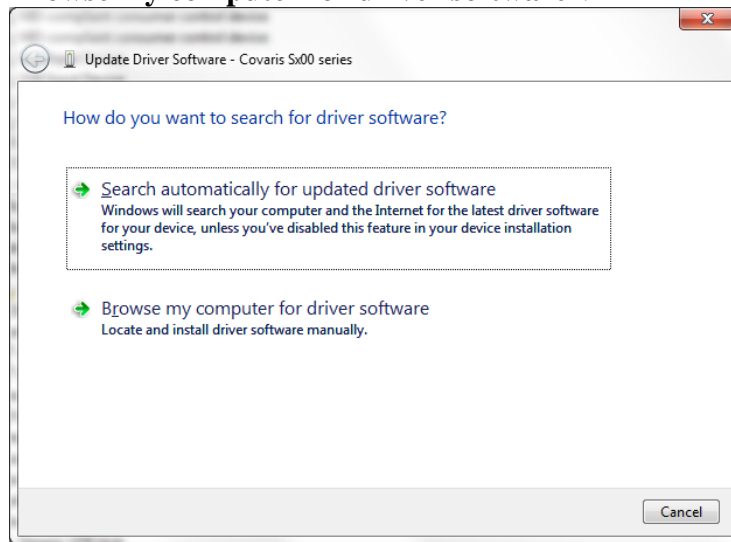


Browse to your desktop and select the **210015 Covaris Driver** folder. Click **Next**.

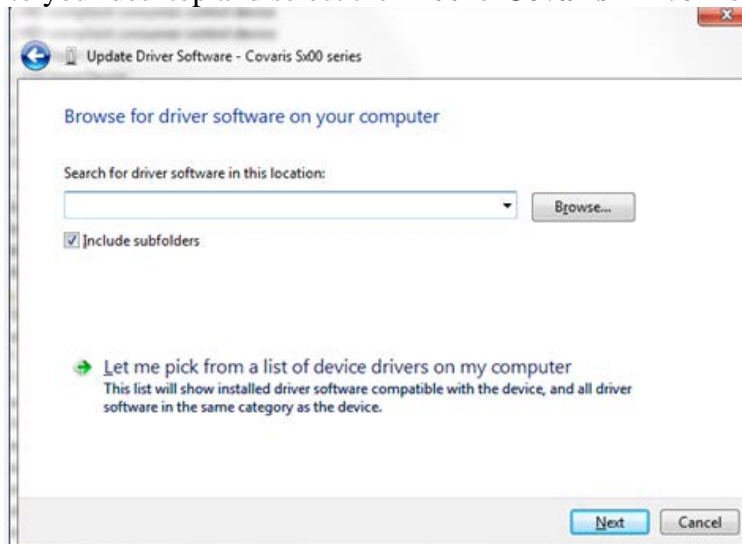

A security warning may appear, as shown below. Click on **“Install this driver software anyway”**

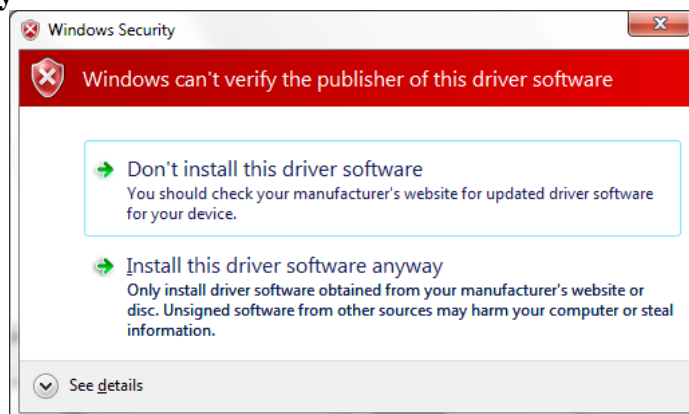


Windows should then install the driver and post the following message. Click on **Close**.

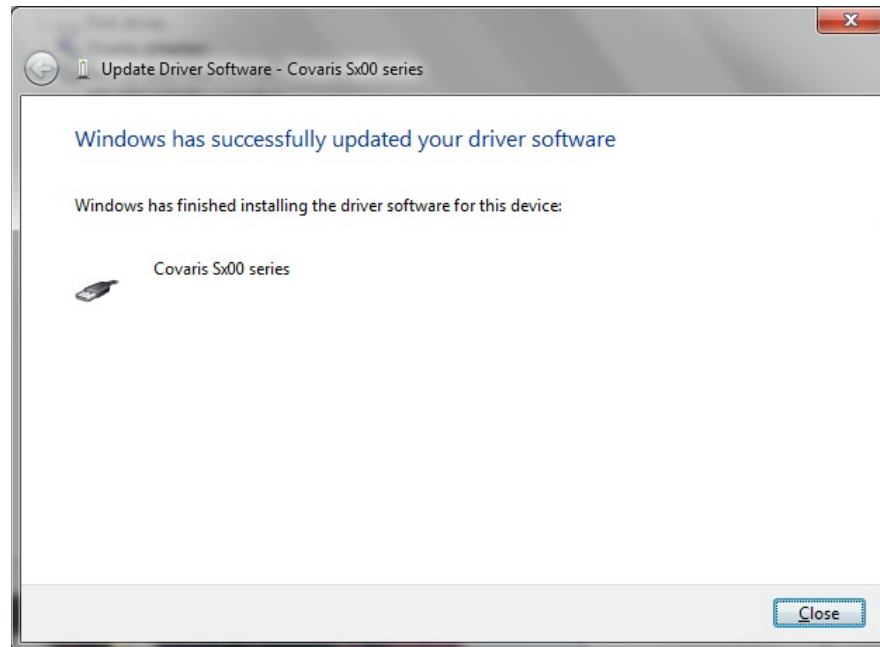

Now return to device manager and under **Other devices**, “USB Serial Port” will appear as below.

Right click on **USB Serial Port** and select **Update driver software**. Repeat previous driver installation process.

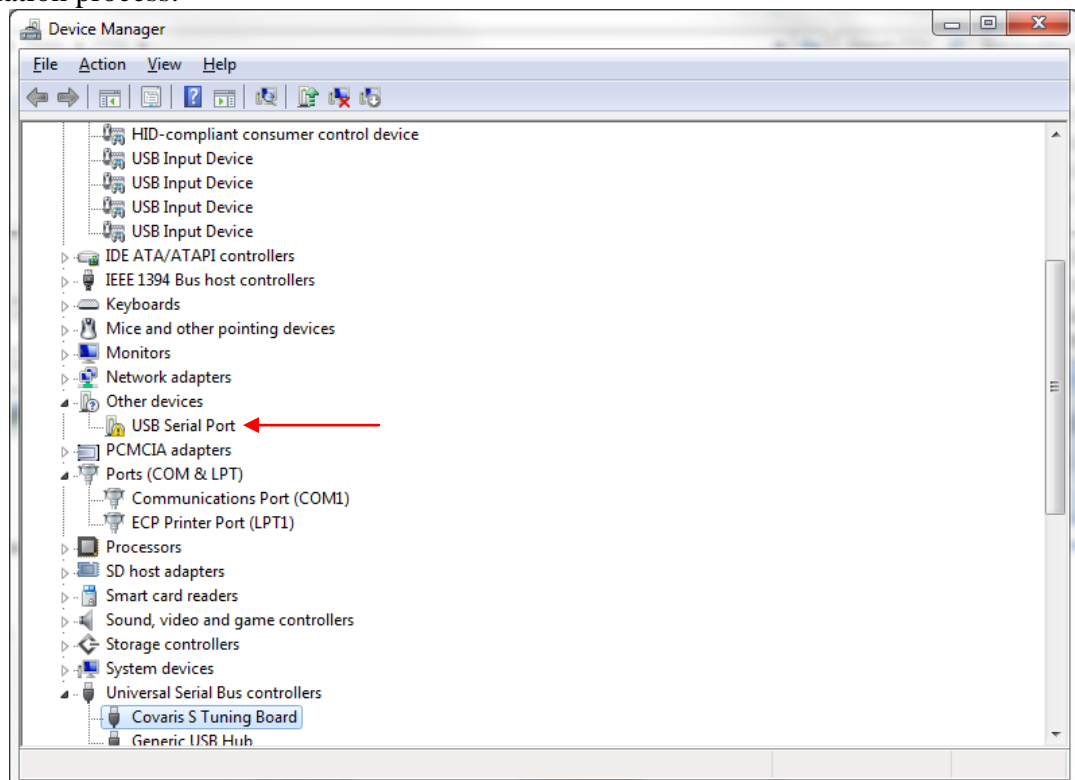

Click **Start Menu**> **Computer**>**C:>Covaris>S-Series vX.X.X>Program**

Right click **SonoLAB v2.x.x** and select **Properties**.

The SonoLAB v2.5.5 Properties window will appear. Click on the **Compatibility tab**.

Under compatibility mode, check off Run this program in compatibility mode for: **Select in the drop down menu Windows XP (Service Pack 2)**

Click **Apply** and **Ok**.

Right click on the SonoLAB S-Series vX.XX and send to **Desktop**.